



BioLab 3.1 Quick Start



BIO LAB 3.1 QUICK START GUIDE

Contact/Customer Support	2
Safety	2
BioNex Software Installation Guide	3

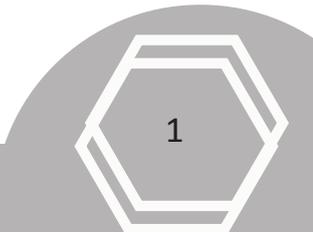

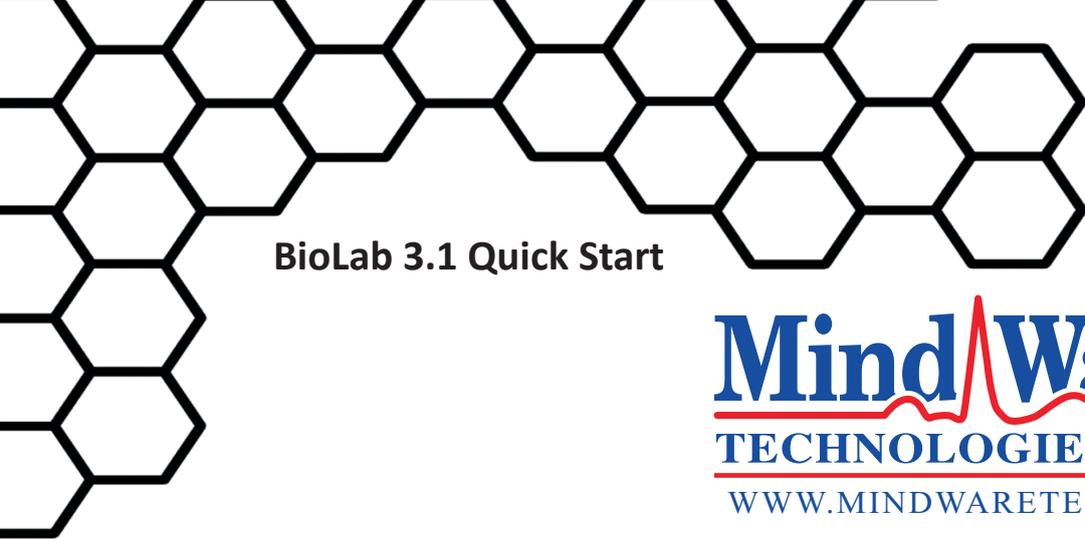

Customer Support:
Email: support@mindwaretech.com
Phone: (614) 626-4888
US Toll Free: (888) 765-9735
Fax: (614) 626-4915

BioLab 3.1 Quick Start



Internet Support

E-mail: info@mindwaretech.com
Web Address: www.mindwaretech.com

Phone


Ph: (614) 626-4888
Fax: (614) 626-4915
MindWare Technologies LTD.
1020-F Taylor Station Rd. • Gahanna, OH 43230

Copyright 2014 © by MindWare Technologies LTD. All Rights Reserved.


Safety

Warning Regarding Medical and Clinical Use of MindWare Technologies LTD Products.

MindWare Technologies LTD products are not designed with components and testing for a level of reliability suitable for use in treatment and diagnosis of humans or as critical components in any life support systems whose failure to perform can reasonably be expected to cause significant injury to a human. Applications of MindWare Technologies LTD products involving medical or clinical treatment can create a potential for death or bodily injury caused by product failure, or by errors on the part of the user or application designer. Any use or application of MindWare Technologies LTD products for or involving medical or clinical treatment must be performed by properly trained and qualified medical personnel, and all traditional medical safeguards, equipment, and procedures that are appropriate in the particular situation to prevent serious injury or death should always continue to be used when MindWare Technologies LTD products are being used. MindWare Technologies LTD products are NOT intended to be a substitute for any form of established process, procedure, or equipment used to monitor or safeguard human health and safety in medical or clinical treatment.

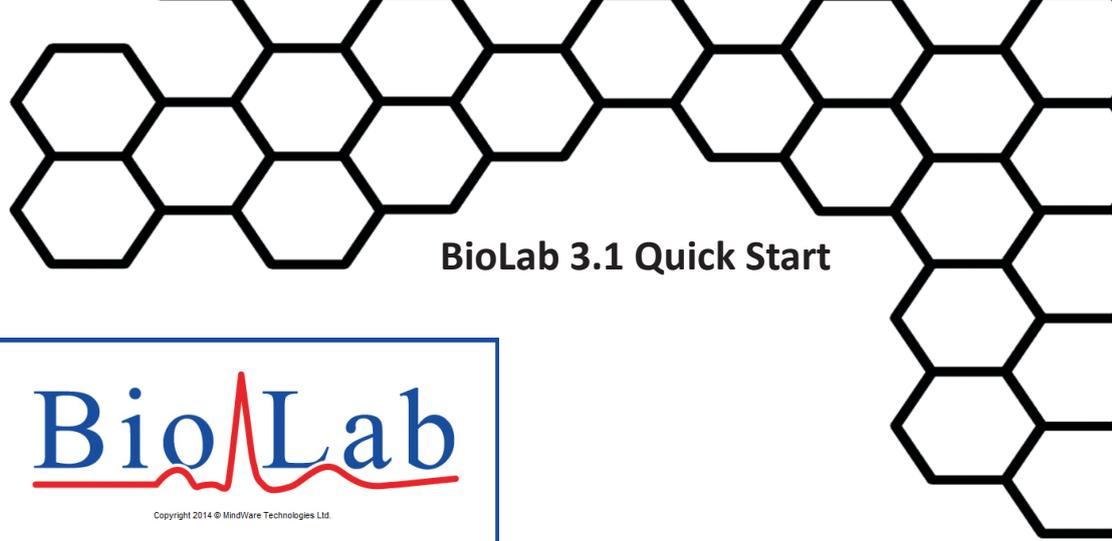

BioLab 3.1 Quick Start


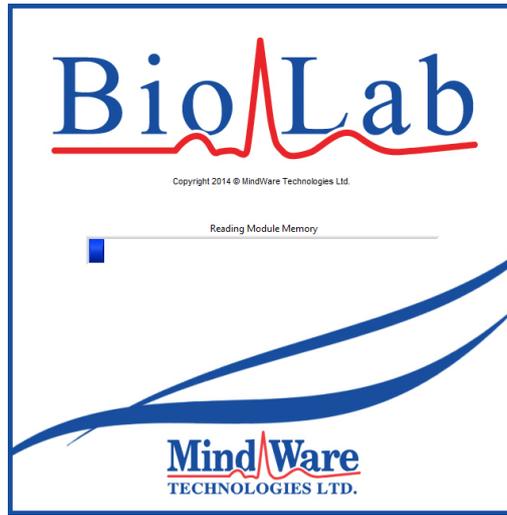


BioNex Software Installation Guide


Note: To ensure proper installation, do not plug the BioNex hardware in until after all software has been installed.

Locate the setup.exe file on the provided USB Drive (or CD/Download) and double click it to begin installation.

Note: The user installing this software must have administrative rights on the machine receiving the installation.

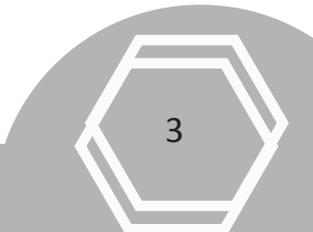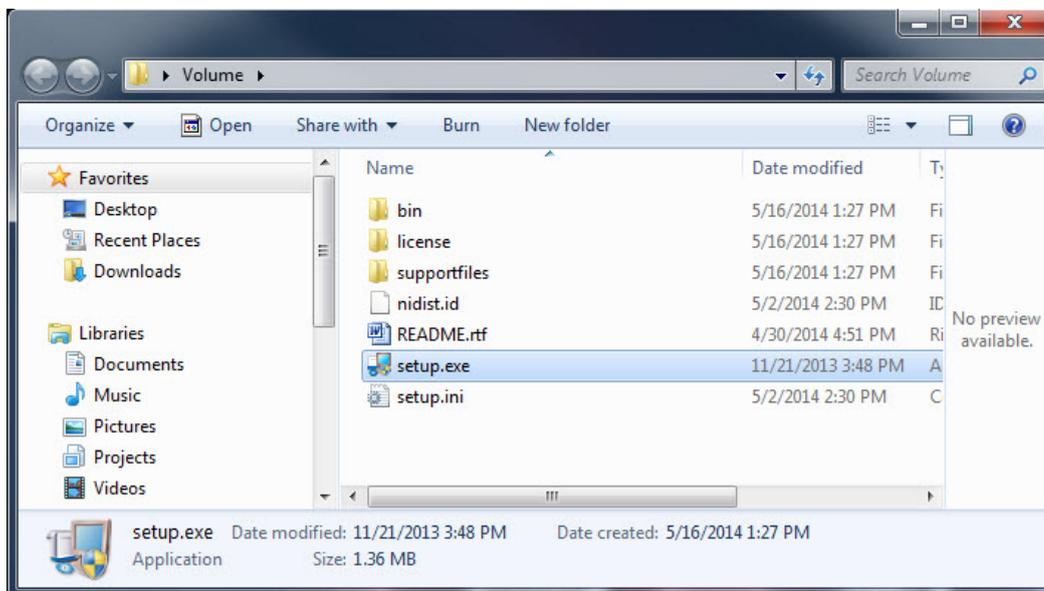


BioLab 3.1 Quick Start

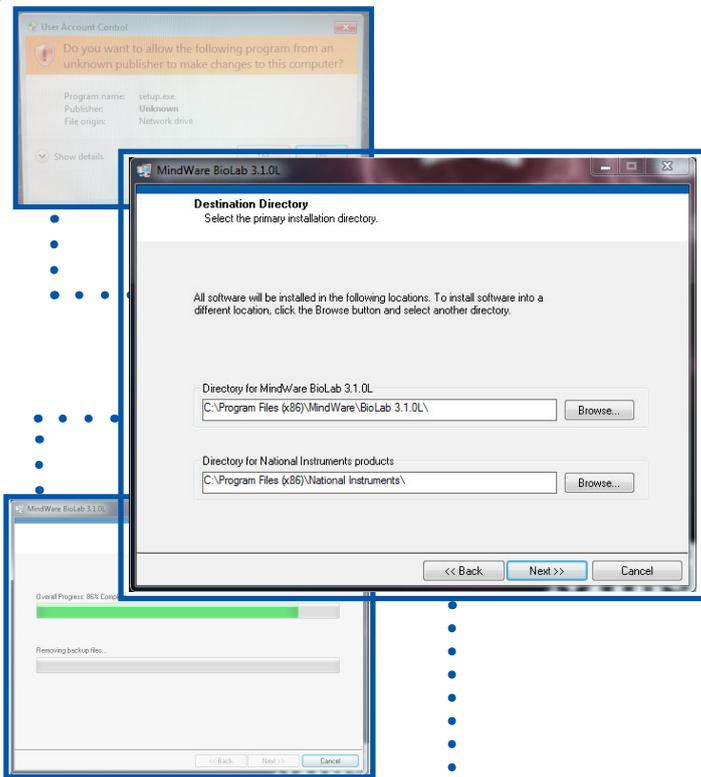

Step 1: Once the installer is run the destination directory will appear indicating the location that the software files will be stored. Hit “Next” twice to continue.


Note: If pop-up appears when running the installer, choose “Yes” to allow install.

Note: Do not change the default directories

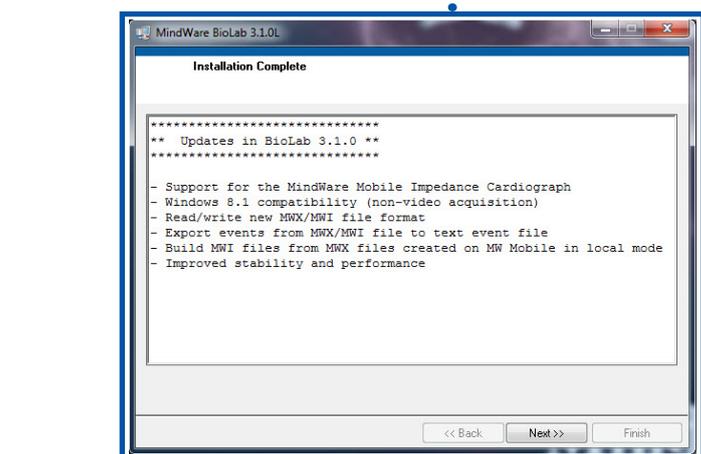


Step 2: Installation Complete screen will appear, click “Next” to continue.

Note: This screen lists the software installed. Please review before continuing.

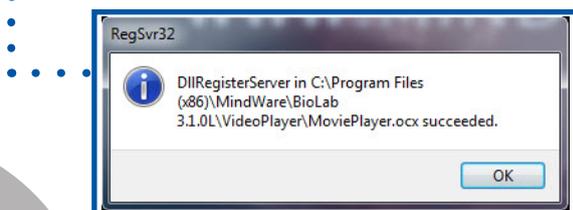


Step 3: Window announcing Video Player installation succeeded will appear, click “OK”.

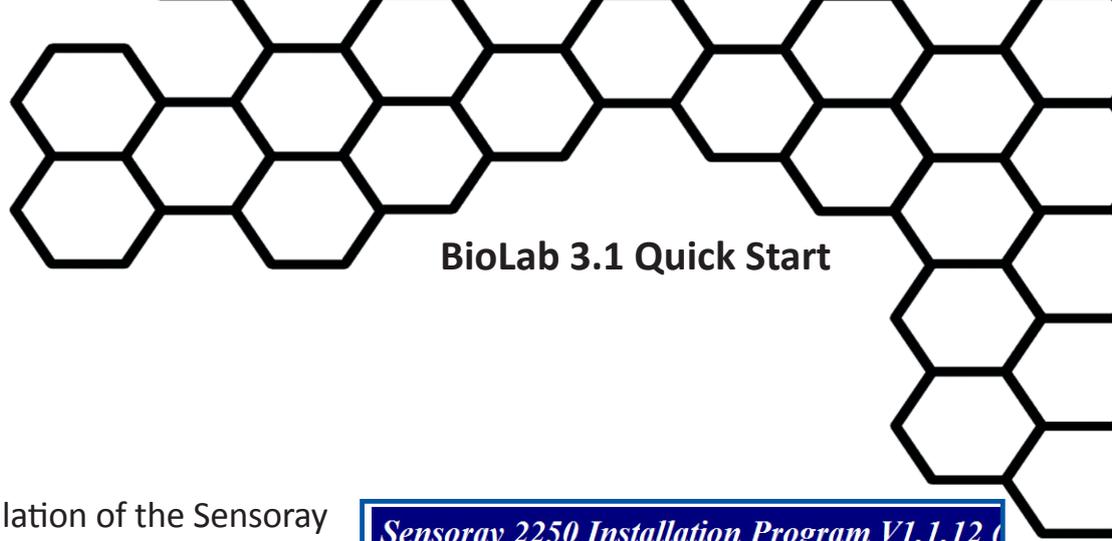

BioLab 3.1 Quick Start

Step 4: This begins the installation of the Sensoray software, click “I Agree” to continue. If you agree you will proceed to the next screen, click “Next” to continue.

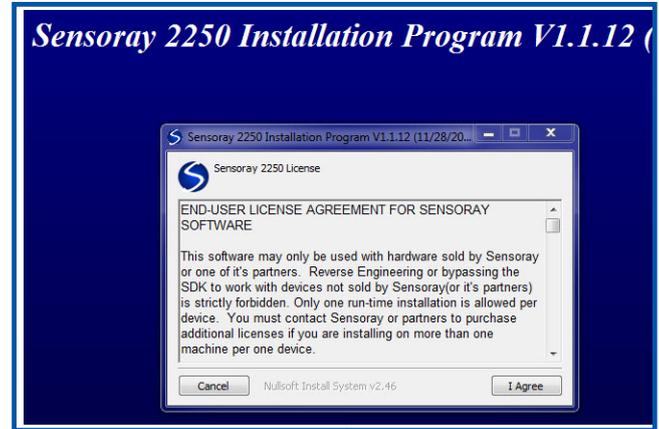





Step 5: This screen shows Sensoray destination directory and begins installation of the program. Click “Install”

*Note: If Sensoray has previously been installed on the computer it will ask that older versions be uninstalled before continuing. Otherwise, continue to **Step 6**.*


Step 5a You may be asked to uninstall, click “Yes” on this screen, then “Uninstall” again on the following screen.

Note: Any prompts that appear asking permission to remove directories should be approved by clicking “Yes”.


Step 5b When uninstall is complete, click the “Close” button in the bottom right of the window.

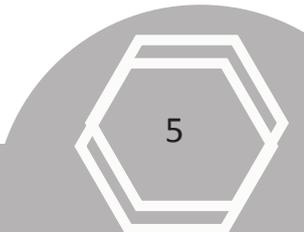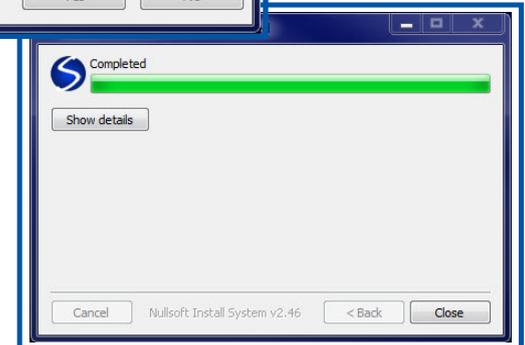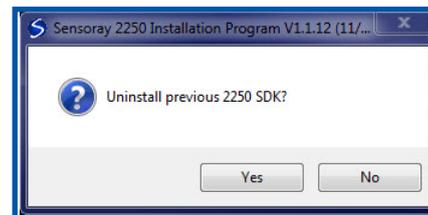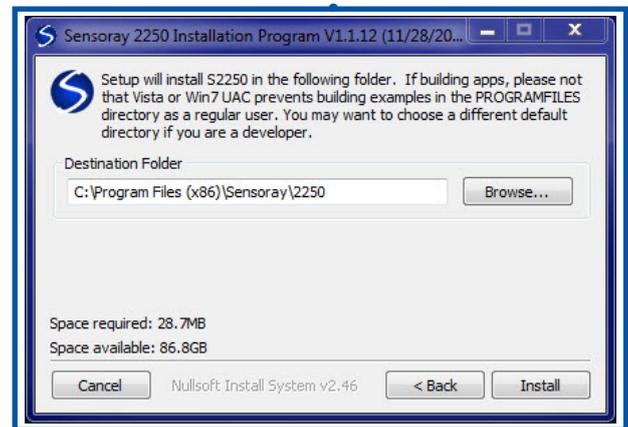


BioLab 3.1 Quick Start

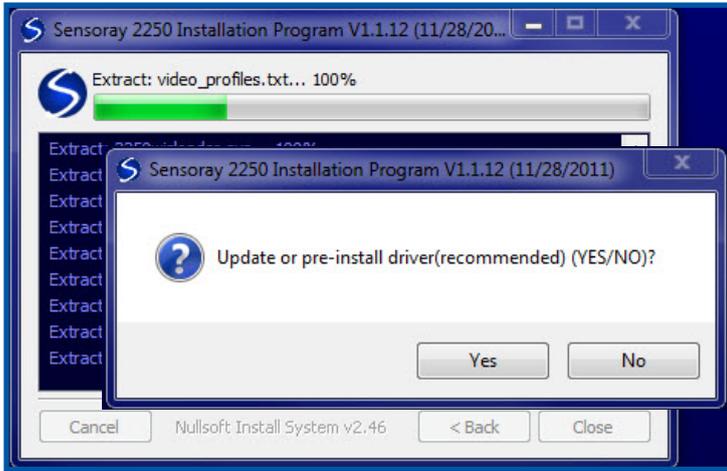


Step 6: Next window will ask you to update your driver , click “Yes” then “OK”.

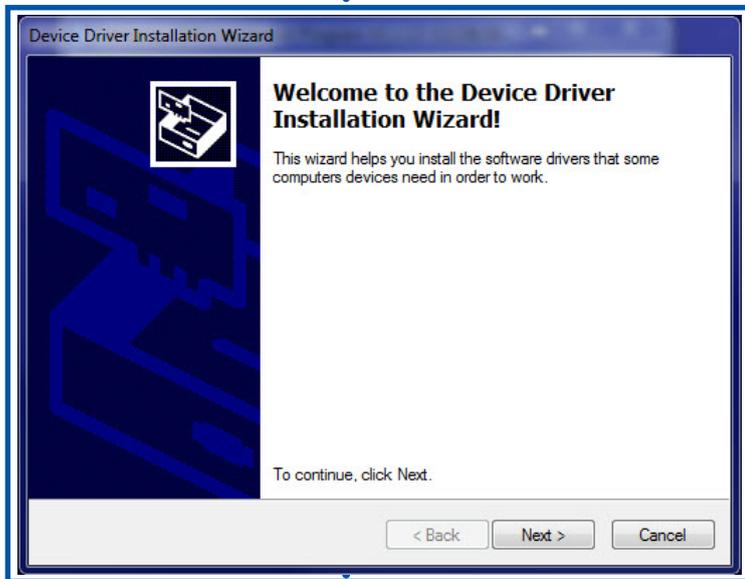

Step 7: Device Driver Installation window will appear. Click “Next” then “Finish”.

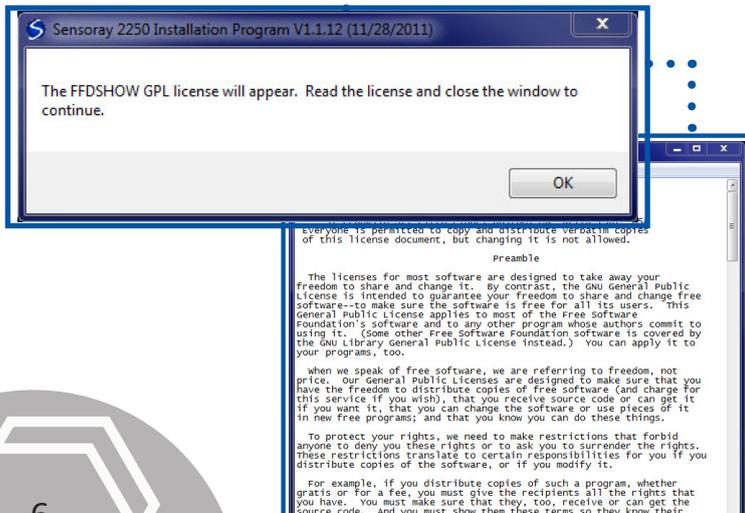

Step 8: Window will prompt you to read the license agreement. Click “OK” then once the agreement has been reviewed, click the “X” in upper right hand corner of the window to close.

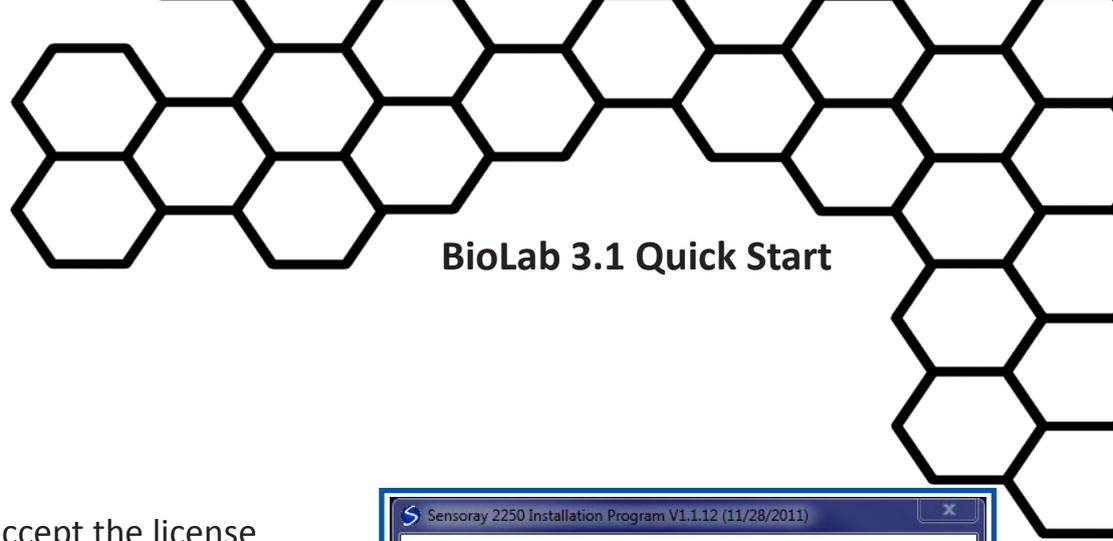


BioLab 3.1 Quick Start

Step 9: A window asking you to accept the license from previous screen will appear. Click “Yes” to continue.

Note: If you click “No” installer will be canceled.

Step 10: Leave default language to English and click “OK”

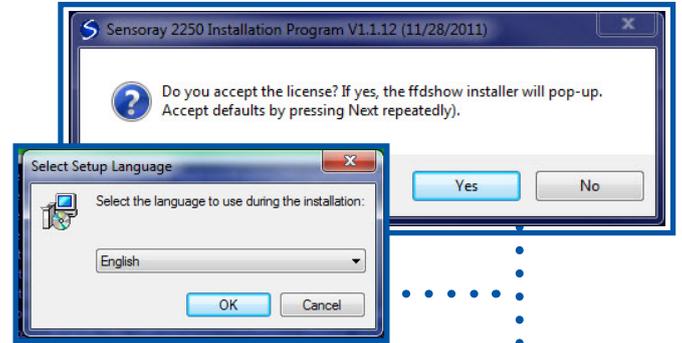

Step 11: On the next few setup screens, accept the defaults by pressing “Next” repeatedly then “Install”.

Note: Leave all settings on default.




Step 12: Completing the setup defaults brings you to this screen. Click “Finish”.

Note: Leave all configuration boxes unselected.



BioLab 3.1 Quick Start

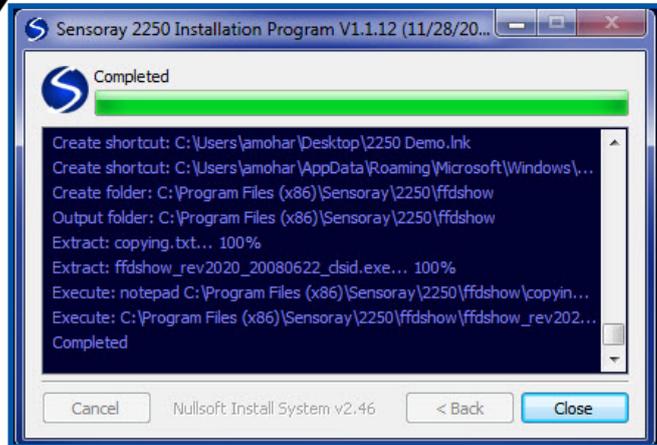

Step 13: Hit “Close” on the completed Sensoray screen.




Step 14: Another Sensoray License Agreement will be shown. To continue click “I Agree”.

Note: If you do not agree the installation will be canceled.


Step 15: In next window click “Next” then “Install”.

Note: Do not change defaults of program file locations.

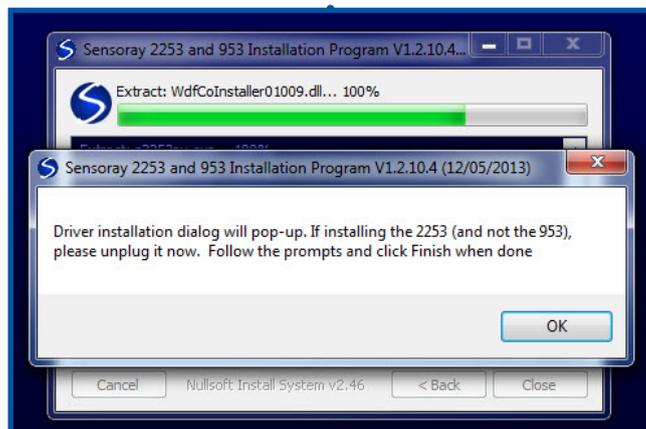


Step 16: On Driver installation dialog pop-up, click “OK”

Note: If instead it asks to Uninstall previous Sensoray driver, click “Yes” then “Uninstall” and “Yes” again to remove directory. Close uninstaller when finished.

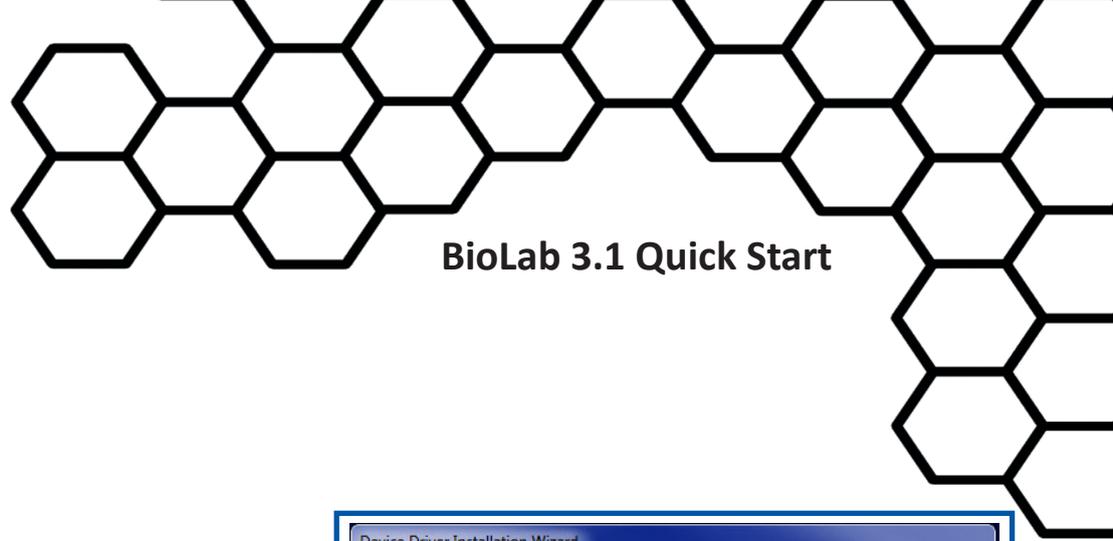

BioLab 3.1 Quick Start


Step 17: Install Wizard will appear. Click “Next” then “Finish”.

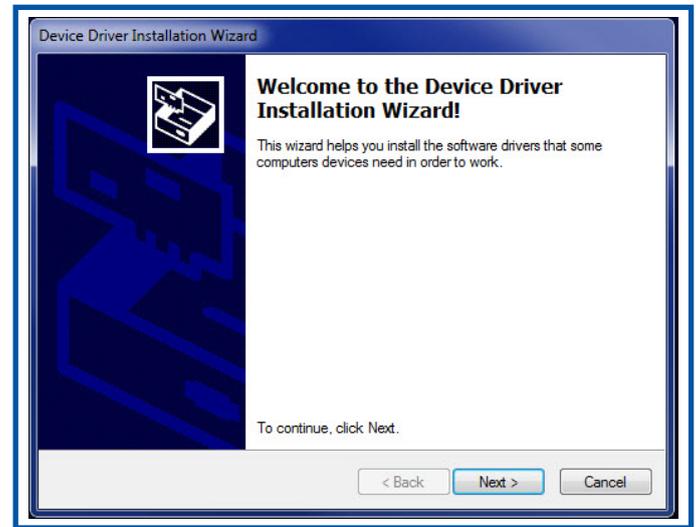

Step 18: Installer will ask wether your are installing 953 or 2253. Choose “No” for 2253.

Note: It is important to select “No” on this screen for installation to complete properly.

Step 19: Click “Close” on installation program window.

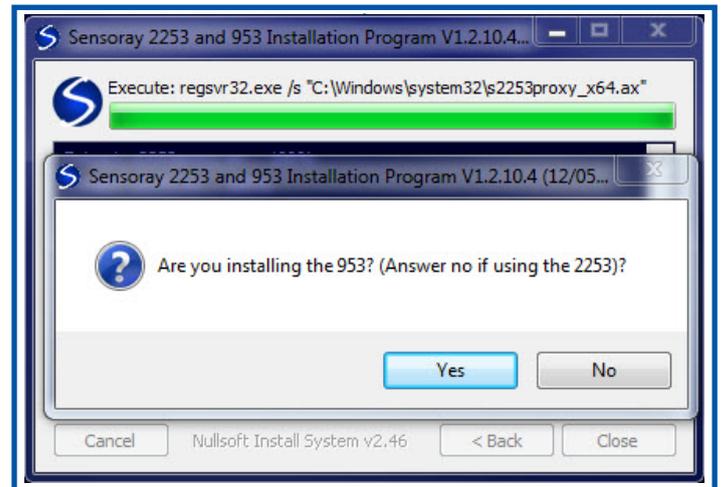

Step 20: Installation is now complete. BioLab 3.1 can now be found in the Start menu.

