

Qualisys Integration Kit

Product Description

Product Description Qualisys Integration Kit

Version 1.0

03/2018

All rights reserved.

Copyright © Tobii AB (publ)

The information contained in this document is proprietary to Tobii AB. Any reproduction in part or whole without prior written authorization by Tobii AB is prohibited.

Products that are referred to in this document may be either trademarks and/or registered trademarks of the respective owners. The publisher and the author make no claim to these trademarks.

While every precaution has been taken in the preparation of this document, the publisher and the author assume no responsibility for errors or omissions, or for damages resulting from the use of information contained in this document or from the use of programs and source code that may accompany it. In no event shall the publisher and the author be liable for any loss of profit or any other commercial damage caused or alleged to have been caused directly or indirectly by this document.

Content subject to change without notice.

Please check Tobii web site www.tobiipro.com for updated versions of this document.

1 Introduction

1.1 Overview

The Qualisys Integration Kit is an extension for the Pro Glasses 2 Head Unit equipped with Qualisys' reflective markers and a connection/synchronization cable which makes it possible to use Pro Glasses 2 together with Qualisys' system.



Please note that all images except the one above shows the kit without the Qualisys reflective markers. Please handle the kit gently to avoid damaging any parts of it.



It is possible to use Tobii Pro Prescription Lenses together with the Qualisys Integration Kit. See Tobii Pro Glasses 2 User's Manual for more information about installing and using prescription lenses.

1.2 Package Content


Item	Quantity
Head unit adapter with reflective markers	1 pc.
Connection/synchronization cable	1 pc.

2 Using the Qualisys Integration Kit

2.1 Mounting the Qualisys Integration Kit onto the Pro Glasses 2 Head Unit

To be able to fit the integration kit onto the head unit you must remove the protective lens mounted on the head unit.


1. Locate the small hook that keeps the lens in place, located between the scene camera and the nose pad at the front of the head unit.
2. Push the hook slightly downwards while gently pulling the lens past the retaining clip.
3. Take the integration kit and place its bottom slot onto the hook above the nose pad and fold the unit upwards into place and finally press lightly on the two top hooks of the unit so it snaps into place.

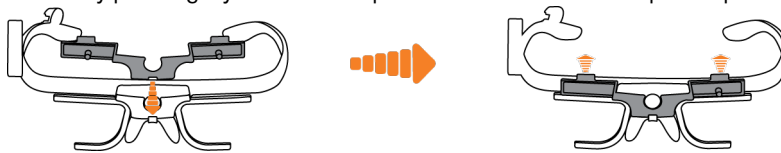

2.2 Removing the Qualisys Integration Kit from the Pro Glasses 2 Head Unit

To remove the integration kit, use the following procedure:

1. Lift the two top hooks off the head unit and gently pull the kit upwards to remove it.

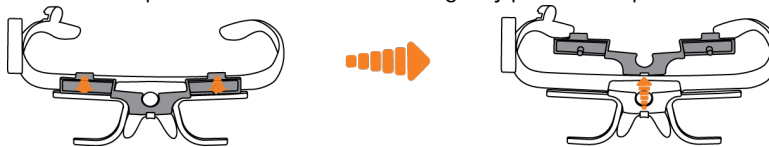

2. Mount the protective lens by hooking the top part of it in the small holes up top and then gently pushing it into place.

2.3 Connecting to Qualisys hardware/software

Use the enclosed connection/synchronization cable for connection to Qualisys' hardware. More information about connecting to and using Qualisys' hardware and software, see the corresponding Qualisys documentation.



©Tobii®. Illustrations and specifications do not necessarily apply to products and services offered in each local market. Technical specifications are subject to change without prior notice. All other trademarks are the property of their respective owners.

Tobii Pro Support

EUROPE / GLOBAL

Phone (SWE): +46 8 522 950 10
Phone (GER): +49 69 24 75 03 4-27
support@tobii.com
Support hours: 8 am - 6 pm
Between July-August: 9 am - 5 pm
(Central European Time, GMT +1)

NORTH AMERICA

Phone: +1 703 738 1320
support.us@tobii.com
Support hours: 8 am - 8 pm
(US Eastern Standard Time, GMT -6)

JAPAN

Phone: +81 3 5793 3316
support.jp@tobii.com
Support hours: 9 am - 5.30 pm
(Japan Standard Time, GMT +9)

CHINA

Phone: +86 180 1558 5168
support.cn@tobii.com